

## Laser Pointer Stand (Order Code STAND)



This light-weight stand is a perfect support for the Laser Pointer available from Vernier (order code LASER) or any other laser pointer that has a 7/16 inch outside diameter. A pair of clips attached to the pivoting bracket securely holds the laser in place. This assembly allows you to accurately point the laser at a photogate.



- To attach the laser pointer, position the laser pointer such that the ON/OFF push touch button is aligned with one of the clips. Push the laser pointer into the clips so that the pointer snaps into place. This arrangement will depress the laser's ON/OFF button, forcing the laser to be on. To turn the laser off, rotate the barrel of the laser so that the touch button is no longer depressed.
- Unscrew the adjustment knob, adjust the height and angle of the bracket to its desired position, then tighten the knob to secure the bracket.



Vernier Science Education  
13979 SW Millikan Way • Beaverton, OR 97005-2886  
Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440  
info@vernier.com • [www.vernier.com](http://www.vernier.com)

Rev. 5/13/2024

